

EasySort™-5 Magnet

Magnet for column-free immunomagnetic separation

Cat.No: EC001

Quantity: 1Unit

Cat.	Products	Holds	Working volume
EC001	EasySort™ Magnet	6+1	0.1-1 mL
	Manual		One Copy

Properties

Storage: Store at ambient temperature of 15-25°C.

STRONG MAGNETIC FIELD: Please keep away from pacemakers, magnets, electronic products such as computers and watches, and other objects that respond to magnetic fields.

Introduction

Elabscience's EasySort™ Magnet is a magnetic separation rack for immune cell sorting, also known as magnetic rack, magnetic sorting rack, etc., which can be used with Elabscience's EasySort™ Cell Isolation Kits.

The EasySort™ Magnet is cylindrical, approximately 65 mm in height and 75 mm in diameter, with a mass of 595 ± 5 g. Each EasySort™ Magnet contains 7 separation holds, each of which can be used for cell separation in one 2 mL round-bottomed microcentrifuge tube or a 5 mL flow tube, and can isolate 0.1-1 mL of cell suspension at a time.

For detailed cell magnetic sorting procedures, please refer to the product manual of EasySort™ Cell Isolation Kits.



Cautions

1. Please keep the product clean and dry, After each time of using, please use pure water to wipe the residual waste liquid or impurities, and use 75% ethanol to scrub the surface to disinfect, then dry it as soon as possible after disinfection.
2. Please avoid contacting this magnet with certain buffers or reagents that have metal corrosive effects to ensure the good performance of this magnet.
3. This magnet has a strong magnetic field, please keep away from pacemakers, computers, cell phones, watches and other electronic products to prevent the dysfunction of related products.

For Research Use Only