

# 2-step Plus Poly-HRP Anti Mouse IgG Detection System (with DAB solution)

Cat. No: E-IR-R216

Size: 3 mL/ 6 mL/ 18 mL



## Product Content

| Cat           | product                                      | 3 mL   | 6 mL   | 18 mL           | Storage |
|---------------|--|--------|--------|-----------------|---------|
| E-IR-R216A    | Normal Goat Blocking Buffer (Ready-to-Use)   | 3 mL   | 6 mL   | 18 mL           | 2~8 °C  |
| E-IR-R216B    | Polyperoxidase-anti-Mouse IgG (Ready-to-Use) | 3 mL   | 6 mL   | 18 mL           | 2~8 °C  |
| E-IR-R216C    | 3% H <sub>2</sub> O <sub>2</sub>             | 3 mL   | 6 mL   | 18 mL           | 2~8 °C  |
| E-IR-R216D    | DAB Concentrate (20×)                        | 150 μL | 300 μL | 900 μL          | 2~8 °C  |
| E-IR-R216E    | DAB Substrate                                | 3 mL   | 6 mL   | 18 mL           | 2~8 °C  |
| <b>Manual</b> |  |        |        | <b>One Copy</b> |         |

## Introduction

2-step plus is a two-step immunohistochemical broad spectrum detection reagent. It polymerizes monovalent Fab fragments of secondary antibody and enzyme, which replaces the secondary antibody and tertiary antibody in traditional method, can directly amplify the binding signal of antibody-antigen. This method not only retains the specific binding ability of antibody with antigen, but also can effectively avoid space steric hindrance caused by excessive polymer molecules. Compared with the traditional SP three-step method, this kit has the characteristics of simple, rapid, high-sensitivity. This system abandons the using of biotin, so it can avoid background staining by endogenous biotin. It can be used in IHC, in which the primary antibody is monoclonal/polyclonal antibody derived from Mouse.

Diluent for DAB concentrated solution has been provided in this kit to avoid the influence of the different water acidity and alkalinity on the DAB Chromogenic Agent.

## Sample dyeing

1. Dewax and hydrate the paraffin section.
2. Make thermal repair or digestion treatment to antigen of the tissue section if necessary according to antigen/antibody situation.
3. Incubate with E-IR-R216C (3% H<sub>2</sub>O<sub>2</sub>) for 10 min to eliminate endogenous peroxidase activity. Wash with PBS or TBS, 2 min×3 times.
4. Add E-IR-R216A (Normal Goat Blocking Buffer (Ready-to-Use)), incubate at 37 °C for 30 min. Shake off any excess liquid.
5. Add primary antibody (Mouse-IgG) with proper dilution ratio, incubate at 20~37 °C for 1~2h or at 4 °C overnight (then rewarm at 37 °C for 30 min). Wash with PBS or TBS, 2 min×3 times. Dry the section with absorbent paper.
6. Add E-IR-R216B (Polyperoxidase-anti-Mouse IgG), incubate at room temperature or 37 °C for 20 min. Wash with PBS or TBS, 2 min×3 times.
7. Add 1 drop (approximately 50 μL) of E-IR-R216D (DAB Concentrate) into each 1 mL of E-IR-R216E (DAB Substrate), mix fully and the mixed reagent is the DAB Working Solution. Prepare fresh solution before use and the prepared solution should be stored in the dark. Fresh prepared DAB Working Solution is valid within 4 hours and the unused solution must be abandoned
8. Take control of the DAB coloration period, the color of tan or brownish yellow is the positive signal. Avoid of excessive reaction.
9. Wash the section with deionized water terminate the chromogenic reaction, then operate the procedures of counterstaining, dehydrating, transparentizing and sealing

## Storage

Store at 2~8 °C, shading light. Avoid of freezing. Valid for 12 months. The reagents are valid within 6 months after opening.

## For Research Use Only

A Reliable Research Partner in Life Science and Medicine

Toll-free: 1-888-852-8623

Web: [www.elabscience.com](http://www.elabscience.com)

Tel: 1-832-243-6086

Email: [techsupport@elabscience.com](mailto:techsupport@elabscience.com)

Fax: 1-832-243-6017

Rev. V1.3