

EasySort™ Human B Cell Isolation Kit

Cat. No: MIH004N

Size: 10 Assays/100 Assays/200 Assays

Component	Component Name	10 Assays	100 Assays	200 Assays	Storage
MIH004NA	EasySort™ Human B Beads Streptavidin 1.0-N	270 µL	1.35 mL×2	1.35 mL×4	2-8°C
MIH004NB	EasySort™ Human B Cell Isolation Cocktail	325 µL	1.65 mL×2	1.65 mL×4	2-8°C
	Manual			1 copy	

Storage

Store at 2-8°C with shading light for 1 year. Avoid freezing and thawing.

Description

Human B cell isolation kit utilizes a negative selection strategy to isolate B cells from fresh or frozen human PBMC sample. The principle of this kit is to use different biotinylated monoclonal antibodies labeling non-B cells, followed by streptavidin conjugated magnetic beads incubation. Unwanted labeled cells are efficiently removed by magnetic isolation, and high purity B cells are isolated.

EasySort™ Human B Cell Isolation Kit can help researcher isolate high purity human B cells with simple experimental procedure. The kit is suitable for isolation of B cells from fresh human PBMC or frozen PBMC, and the isolated B cells can be directly used for downstream applications. The B cells isolated from normal PBMC using this kit is typically $94.1 \pm 2.6\%$.

Reagents and Materials Not Supplied

1. Reagents:

PBS, fetal bovine serum (FBS), EDTA, Human peripheral blood mononuclear cells separation solution, DNase I

2. Materials:

70 µm mesh nylon strainer, 1.5 mL/2 mL EP tube, 15 mL/50 mL centrifuge tube, flow tube

3. Instrument:

Optical microscope, centrifuge, 5 mL magnetic rack

For Research Use Only

Experimental Operation

NOTE: The following operations must be performed under sterile conditions

➤ Isolation buffer preparation

Add fetal bovine serum (final concentration of 2%) and EDTA (final concentration of 2 mM) to PBS buffer and filter the prepared buffer with 0.22 µm filter.

NOTE: Sealed store the prepared buffer at 4°C and use within 1 week. In addition, 2% fetal bovine serum can be replaced by 0.5% BSA.

➤ Sample Preparation and Processing

1. Fresh human PBMC: PBMC sample is obtained from fresh human whole blood by density gradient centrifugation. Wash PBMC twice with isolation buffer, centrifuge at 300 g for 5 min, filter the PBMC through a 70 µm mesh nylon strainer and adjust the cell density to 1×10^8 cells/mL for cell isolation.

Note: The best separation effect can be achieved when the freshly collected human blood is separated within 1 hour. Approximately 1×10^7 PBMC can be obtained from 10 mL of human blood.

2. Frozen PBMC: incubate the frozen PBMC should be incubated with DNase I solution (PBS) at a concentration of 100 µg/mL for 15 min at room temperature before cell isolation. Wash sample twice with isolation buffer, centrifuged at 300 g for 5 min. Filter aggregated suspensions through a 70 µm mesh nylon strainer and adjust cell density at 1×10^8 cells/mL.

➤ Cell Isolation

- a) Prepare 100 µL of cell suspension (about 1×10^7 cells), add 32.5 µL Human B Cell Isolation Cocktail, mix fully and incubate for 5 min at room temperature.

Note: Please make sure the cells are single-cell suspension.

- b) Add isolation buffer to a final volume of 2 mL, centrifuge at 300 g for 5 min. Discard the supernatant, and then resuspend the cells with 100 µL isolation buffer.
- c) Wash Beads Streptavidin 1.0-N: Vortex beads for 20 seconds, add 27 µL Beads in 1.5 mL EP tube. Put the tube on a 5 mL magnetic rack (self-provided) and stand for 30 seconds. Remove the supernatant, then resuspend beads with 1 mL isolation buffer, and stand for 5 minutes at room temperature. Remove the supernatant, then resuspend beads with 27 µL isolation buffer.
- d) Transfer the cells to the bottom of the flow tube (**Note: Avoid adding along tube walls**), add 27 µL washed Human B Beads Streptavidin 1.0-N, mix gently and incubate at room temperature for 5 min.

Note:

- ✧ If more than 1×10^7 cells are to be isolated, increase the amount of Human B Cell Isolation Cocktail and Human B Beads Streptavidin 1.0-N proportionally while ensuring the cell density remains 1×10^8 cells/mL. If fewer than 1×10^7 cells are to be isolated, resuspend the cells with 100 µL isolation buffer, add 32.5 µL Human B Cell Isolation Cocktail and 27 µL washed Human B Beads Streptavidin 1.0-N.

For Research Use Only


✧ The 5 mL flow tube is suitable for less than 1×10^8 cells.

- e) Add isolation buffer to a final volume of 2.5 mL, mix gently with a pipette by blowing up and down for 7-8 times until no particles of magnetic beads are visible. Put the tube on a 5 mL magnetic rack (self-provided) and stand for 5 min.

Note: Please mix the liquid thoroughly to avoid the magnetic beads clumping and affecting the isolation efficiency.

- f) Transfer the cell suspension to a clean centrifuge tube, centrifuge at 300 g for 5 min. Discard the supernatant, resuspend the cells with buffer required for the subsequent experiments.

Typical data

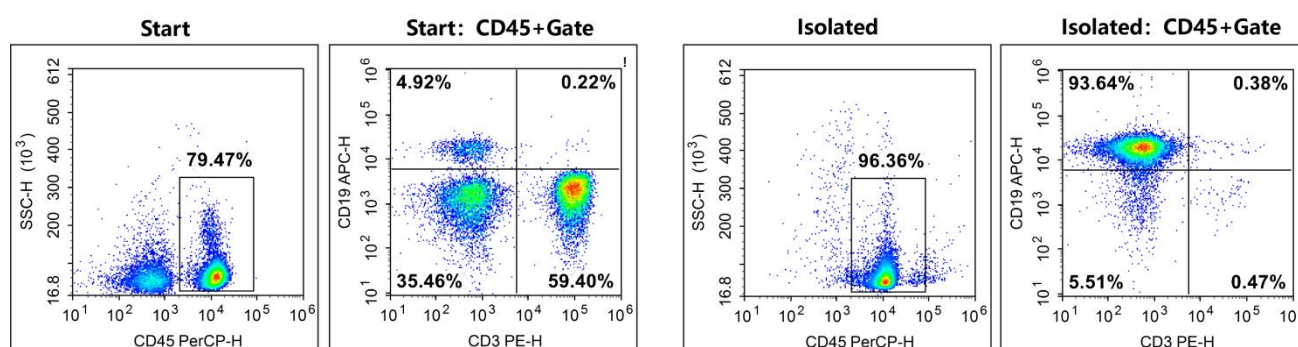

As shown in the figure above, flow cytometric analysis of B cells purity before and after sorting revealed proportions of 4.9% and 93.6%, respectively.

Fluorochrome-conjugated antibody	Cat.	Company
PerCP Anti-Human CD45 Antibody[HI30]	E-AB-F1137F	Elabscience
PE Anti-Human CD3 Antibody[OKT-3]	E-AB-F1001D	Elabscience
APC Anti-Human CD19 Antibody[CB19]	E-AB-F1004E	Elabscience

Cautions

1. This kit is for research use only.
2. Please take safety precautions and follow the procedures of laboratory reagent operation.
3. Avoid freezing and thawing during the use and storage of the beads.
4. The data presented for this product are based on testing of peripheral blood mononuclear cells (PBMCs) from healthy individuals. When processing peripheral blood from non-healthy individuals, to ensure optimal experimental outcomes, first assess the B cells content in the sample. If the B cells content is below 3%, it is recommended to perform a pre-sorting test, appropriately increase the amount of antibodies and beads per assay, or conduct a secondary sorting step.
5. The cell clusters in the cell suspension will affect the purity of cell isolation. Therefore, cell

For Research Use Only

suspension should be filtered with a 70 µm mesh nylon sieve before formal isolation.

6. Cell suspension should be isolated immediately after preparation, the longer the storage time, the greater the impact on cell activity.
7. The cell suspension and reagents should be added directly to the bottom of flow tube to avoid sticking to the wall, resulting in insufficient reaction and affecting the isolation efficiency.
8. In order to ensure the activity of the cells, the whole process of the experiment should be completed on ice as much as possible, except for the incubation at room temperature.
9. It is recommended to use low adsorption pipette tips and centrifuge tubes to avoid the loss of magnetic beads and antibodies due to adsorption.
10. The kit should be used in combination with a magnetic rack.